



CELL LAB Rat Anti-Mouse IgE (ϵ chain specific)

<u>Cat. No.</u>	<u>Form</u>	<u>Quantity</u>
732392	Purified (UNLB) Antibody	0.5 mg
732393	Fluorescein (FITC) Conjugate	0.5 mg
732394	Alkaline Phosphatase (AP) Conjugate	1.0 mL
732395	Horseradish Peroxidase (HRP) Conjugate	1.0 mL
732396	Biotin (BIOT) Conjugate	0.5 mg
732397	Phycoerythrin (PE) Conjugate	0.1 mg
733236	Phycoerythrin (PE) Conjugate	0.2 mg

For Laboratory Use Only

DESCRIPTION

Clone: 23G3
Isotype: Rat IgG1 κ
Specificity: Reacts with the ϵ heavy chain of BALB/c mouse IgE as demonstrated by ELISA.

APPLICATIONS

- Identification and enumeration of IgE⁺ cells by flow cytometry
- Identification and enumeration of IgE⁺ cells by immunofluorescence microscopy
- Enzyme-Linked Immunosorbent Assay (ELISA)

CHARACTERIZATION

To ensure lot-to-lot consistency, each batch of product is tested to conform with characteristics of a standard reference reagent using ELISA, Particle Concentration Fluorescence Immunoassay and/or flow cytometry.

WORKING DILUTIONS

Flow Cytometry:

FITC conjugate	$\leq 1 \mu\text{g}/10^6$ cells
BIOT conjugate	$\leq 1 \mu\text{g}/10^6$ cells
PE conjugate	$\leq 0.2 \mu\text{g}/10^6$ cells

ELISA:

AP conjugate	1:500-1:1,000
HRP conjugate	1:1,000-1:2,000
BIOT conjugate	1:5,000-1:10,000

Other Applications: Since applications vary, determine the optimum working dilution of the product that is appropriate for your specific needs.

HANDLING AND STORAGE

- The purified (UNLB) antibody is supplied as 0.5 mg of purified immunoglobulin in 1.0 mL of 100 mM borate buffered saline, pH 8.0. No preservatives or amine-containing buffer salts added.
- The fluorescein (FITC) conjugate is supplied as 0.5 mg in 1.0 mL of PBS/NaN₃.

- The alkaline phosphatase (AP) conjugate is supplied as 1.0 mL of stock solution in 50 mM Tris/1 mM MgCl₂/50% glycerol, pH 8.0, containing 0.1% NaN₃ as preservative.
- The horseradish peroxidase (HRP) conjugate is supplied as 1.0 mL of stock solution in 50% glycerol/50% PBS, pH 7.4. No preservative added.
- The biotin (BIOT) conjugate is supplied as 0.5 mg in 1.0 mL of PBS/NaN₃.
- The phycoerythrin (PE) conjugates are supplied as 0.1 mg in 1.0 mL or 0.2 mg in 2.0 mL of PBS/NaN₃ and a stabilizing agent.
- Protect fluorochrome-conjugated forms from light. Do not freeze.
- Reagent is stable until the expiration date on the vial when stored at 2-8°C.

STATEMENT OF WARNINGS

1. Specimens, samples and all material coming in contact with them should be handled as if capable of transmitting infection and disposed of with proper precautions.
2. Never pipet by mouth and avoid contact of samples with skin and mucous membranes.
3. Do not use reagent beyond the expiration date on the vial label.
4. Minimize exposure of reagent to light during storage or incubation.
5. Avoid microbial contamination of reagent or erroneous results may occur.
6. Use Good Laboratory Practice (GLP) when handling this reagent.
7. Harmful if swallowed.
8. After contact with skin, wash immediately with plenty of water.
9. Contains sodium azide. Sodium azide under acidic conditions yields hydrazoic acid, an extremely toxic compound. Azide compounds should be flushed with running water while being discarded. These precautions are recommended to avoid deposits in metal piping in which explosive conditions can develop. If skin or eye contact occurs, immediately wash excessively with water.

TRADEMARKS

The Beckman Coulter logo is a trademark of Beckman Coulter, Inc.

For additional information or if damaged product is received, contact your local Beckman Coulter Representative.



Manufactured for:
Beckman Coulter, Inc.
4300 N. Harbor Blvd.
Fullerton, CA 92835
www.beckmancoulter.com

Printed in USA
Made in USA

© 2005 Beckman Coulter, Inc.
All Rights Reserved.

PN 733867-B